

## **Product Data Sheet**

Upright 16 cu. ft. Glass Door Refrigerator, High Performance - Certified to NSF/ANSI 456 Standard for Vaccine Storage

# PH-ABT-NSF-S16G

#### **Product Description**

These cutting-edge pharmacy refrigerators are certified in accordance with the NSF/ANSI 456 Standard for Vaccine Storage. With this certification, units protect pharmaceuticals at optimal temperatures, preventing waste and allowing for peak delivery. Our Standard line provides multi-functional features in a cost-effective design.

These glass door refrigerators utilize microprocessor controllers and feature temperature alarms, remote alarm contacts, and probe access ports with included probes. Units run on natural, hydrocarbon refrigerant for environmental health and energy efficiency.

General Description and Application				
Description	Single Glass Door Pharmacy/Vaccine Upright Refrigerator			
Operational environment	Indoor use only, +18°C to +26°C (+65°F to +78°F), <70% RH			
Storage capacity	16 cu. ft. gross volume			
Door	One swing glass door, self-closing, right hinged, non-reversible, magnetic sealed gasket, ke lock			
Shelves	Four shelves (three adjustable/one fixed) with guard rail on back			
Mounting	3 1/2" Swivel Casters(two locking)			
Interior lighting	Shielded, switched LED lighting, full coverage, balanced spectrum			
Airflow management	Forced Air technology, patent pending			
External probe access	Rear wall port (3/4") dia.			
Insulation	Cabinet is foamed-in-place with EPA compliant high density urethane foam			
Exterior materials	White powder coated steel			
Access control	Pyxis®, Omnicell® and AcuDose RX® compatible			
General warranty	One (1) year parts and labor warranty, excluding display probe calibration			
Compressor warranty	Five( 5) years compressor warranty			
Product Weight	249			
Shipping Weight	289			
Rated Amperage	3			
Power Plug/Power Cord	NEMA 5-15 plug, 8 to 10 ft typical, conforms to UL471 requirements, Vaccine Storage power			
Facility Electrical Requirement	110-120V AC: 15 A (minimum)			
Agency Listing and Certification	Certified with the temperature performance requirements as defined in the NSF/ANSI 456 Standard for Vaccine Storage for all testing scenarios. UL, C-UL, ETL, C-ETL listed and certified to UL471 standard, hydrocarbon refrigerant safety. Energy Star Certified			
Included Accessories	Temperature monitor device (TMD) complies with the current CDC guidelines, with 3 years certification of calibration, "buffered" probe in the product simulated solution, min/max memory, field installable, and visual & audible temp alarm			

# Pharmacy refrigerator/freezer toolkit and temperature logs

## **Refrigeration System**

Compressor Refrigerant Condenser Evaporator Defrost Hermetic, high performance EPA SNAP compliant, R290, propane Fin and tube design, high efficiency fan Fin and tube design, high efficiency fan Cycle optimized, zero energy

Performance

Uniformity<sup>1</sup> (Cabinet air) +/- 0.9°C Stability<sup>2</sup> (Cabinet air) +/- 1.0°C Maximum temperature variation +/-1.2°C (Cabinet air) Temperature rise after 8 sec door Temperature did not exceed 6.5°C at any probe for all required NSF/ANSI 456 testing protocols<sup>3</sup> openings Recovery after 3 min door opening All probes recover to under 8°C within 4.3 min. 1.25 KWh/day<sup>4</sup> Energy consumption 1.97 KWh/day (280BTU/h)4 Average heat rejection Noise pressure level (dBA) 48 or less installed Pull down time to 4°C nominal 30 min operating temp

Controller, Configuration, Alarms and Monitoring							
Controller technology	Parametric, microprocessor, LED display with 0.1°C resolution						
Temperature setpoint range	1°C to 10°C (Controller settings must remain unaltered to ensure thermal performance compliant with NSF/ANSI 456 Standard for Vaccine Storage requirements)						
Display probe	Calibrated, stainless steel						
External alarm connection	State switching remote alarm contacts						
	Visual and audible indicators						
Alarms	High / Low temperature, compliant with alarm requirements defined in the NSF/ANSI 456						
	Standard for Vaccine Storage						
Simulator ballast	Glass bead thermal media						

Performance data acquired at 22°C ambient, using NSF/ANSI 456 compliant validation ballast probes, empty chamber, during stabilized steady state operation and a DAQ sampling rate of one measurement every 10 seconds

1 - Uniformity is defined as the maximum variance in temperature across all probes at any point in time over the testing period


2 - Stability is defined as the maximum variance in temperature experienced by any single probe over the testing period


3 - Temperature performance for all loaded and unloaded door opening protocols, all alarm, controller and probe requirements as defined in the NSF/ANSI 456 standard for vaccine storage


4 - Data per Energy Star test results or equivalent testing and calculation. Heat rejection based on daily averages, not continuous operation. Performance exceeds Energy Star requirements.





\*-one or more of these certifications may apply to this unit.

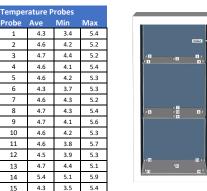

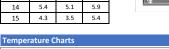

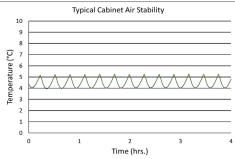

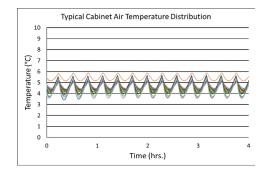

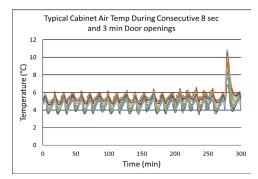



# **Product Data Sheet**


Upright 16 cu. ft. Glass Door Refrigerator, High Performance - Certified to NSF/ANSI 456 Standard for Vaccine Storage







Dimensions										
		Width	Depth	Height	Door Swing	Total open Depth				
	Exterior	25"	26"	79"	23 1/8"	49 1/4"				
	Interior	20"	19 1/8"	48 1/8"						

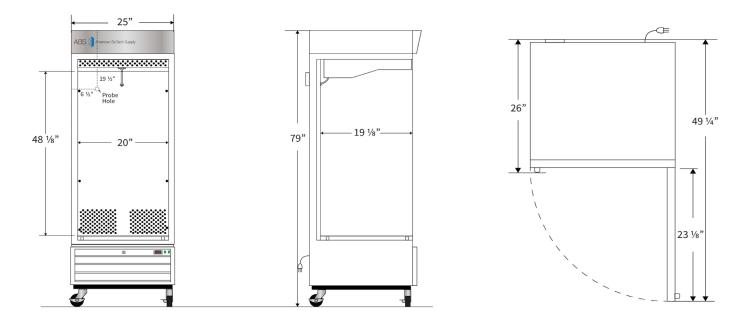

Contact Customer Service Technical Service Rev\_09302022

800-648-4041 Option 3 800-648-4041 Option 5, Parts Option 4 customerservice@horizonscientific.com technicalservice@horizonscientific.com