

PH-ABT-NSF-UCBI-0404G-ADA

Product Description

These cutting-edge pharmacy refrigerators are certified in accordance with the NSF/ANSI 456 Standard for Vaccine Storage. With this certification, units protect pharmaceuticals at optimal temperatures, preventing waste and allowing for peak delivery. Our Premier line includes premium features such as extensive alarm systems and digital touch pad displays.

The ADA compliant, glass door refrigerators utilize microprocessor controllers and feature temperature alarms, remote alarm contacts, LED interior lighting, and probe access ports with included probes. American Biotech Supply Vaccine Storage Refrigerators utilize HFC-free refrigerant for environmental health and energy efficiency.


General Description and Application			
Description	Single Glass Door Pharmacy/Vaccine Undercounter Refrigerator Built-In ADA Compliant		
Operational environment	Indoor use only, +18°C to +26°C (+65°F to +78°F), <70% RH		
Storage capacity	4.6 cu. ft. gross volume		
Door	One swing glass door, self-closing, right hinged, non-reversible, magnetic sealed gasket, keyed lock		
Shelves	Three shelves (two adjustable/one fixed) with guard rail on back		
Mounting	Low profile roller wheels and leveling legs		
Interior lighting	Shielded, switched LED lighting, full coverage, balanced spectrum		
Airflow management	Forced Air technology, patent pending		
External probe access	Rear wall port (1/2") dia.		
Insulation	Cabinet is foamed-in-place with EPA compliant high density urethane foam		
Exterior materials	White powder coated steel		
Access control	Pyxis [®] , Omnicell [®] and AcuDose RX [®] compatible		
General warranty	Two (2) years parts and labor warranty, excluding display probe calibration		
Compressor warranty	Five (5) years compressor warranty		
Product Weight	100 lbs.		
Shipping Weight	140 lbs.		
Rated Amperage	1.74 Amps		
Power Plug/Power Cord	NEMA 5-15 plug, 8 to 10 ft typical, conforms to UL471 requirements, Vaccine storage power cord warning label		
Facility Electrical Requirement	110-120V AC: 15 A (minimum)		
Agency Listing and Certification	Certified in accordance with the NSF/ANSI 456 Standard for Vaccine Storage. UL, C-UL, ETL, C- ETL listed (either single or dual agency listings) and certified to UL471 standard, hydrocarbon refrigerant safety. Energy Star Certified		
Included Accessories	Temperature monitor device (TMD) complies with the current CDC guidelines, with 3 years certification of calibration, "buffered" probe in the product simulated solution, min/max memory, field installable, and visual & audible temp alarm		
	Pharmacy refrigerator/freezer toolkit and temperature logs		

Certifications




*-one or more of these certifications may apply to this unit.

Temperature Probes				
Probe	e Ave Min		Max	
1	4.6	3.5	5.8	
2	4.9	4.3	5.4	
3	5.0	4.4	5.6	
4	4.6	3.4	5.8	
5	5.0	4.6	5.3	
6	5.3	4.7	5.9	
7	4.8	4.2	5.5	
8	5.1	4.5	5.8	
9	4.8	3.9	5.8	
10	4.8	3.9	5.8	
11	5.5	4.9	6.2	
12	5.1	4.6	5.6	
13	4.9	4.3	5.5	
14	4.9	4.0	5.9	
15	5.5	4.9	6.2	

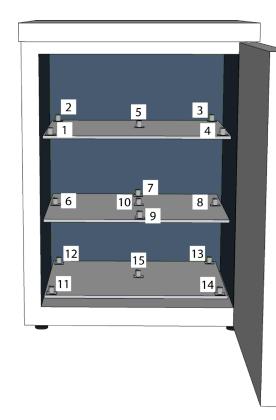

Temperature Charts

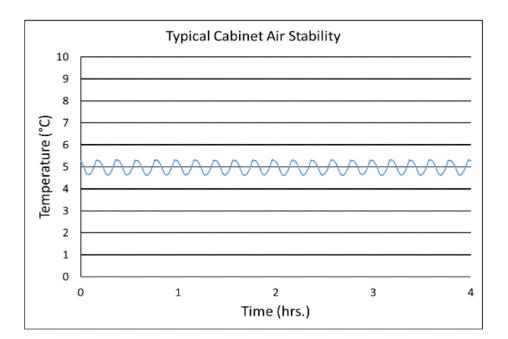

Refrigeration System Compressor Befrigerant EBA SNAB compliant Before and the second secon


Product Data Sheet


Undercounter 4.6 cu. ft. Built-in Glass Door Vaccine Refrigerator ADA - Certified to NSF/ANSI 456 Standard for Vaccine Storage

Reingerant	EPA SNAP compliant, R600a, Isobutane
Condenser	Hybrid fin and tube with low noise fan
Evaporator	Plate wall
Defrost	Cycle optimized, zero energy

Per	forr	nai	nce		
			4	 	

Performance	
Uniformity ¹ (Cabinet air)	+/- 0.8°C
Stability ² (Cabinet air)	+/- 1.2°C
Maximum temperature variation (Cabinet air)	+/- 1.4°C
Temperature rise after 8 sec door openings	Temperature did not exceed 6.4°C at any probe for all required NSF/ANSI 456 testing protocols ³
Recovery after 3 min door opening	All probes recover to under 8°C within 4.8 min.
Energy consumption	1.15 KWh/day⁴
Average heat rejection	1.57 KWh/day (224 BTU/h) ^₄
Noise pressure level (dBA)	43 or less installed
Pull down time to nominal operating	35 min
temp	

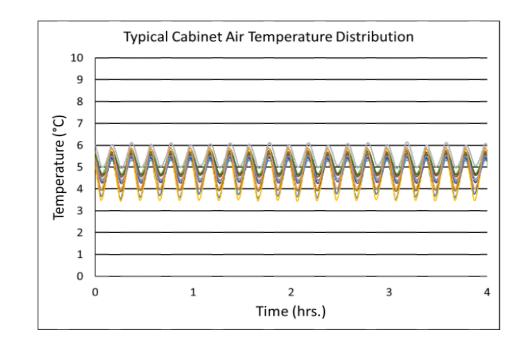

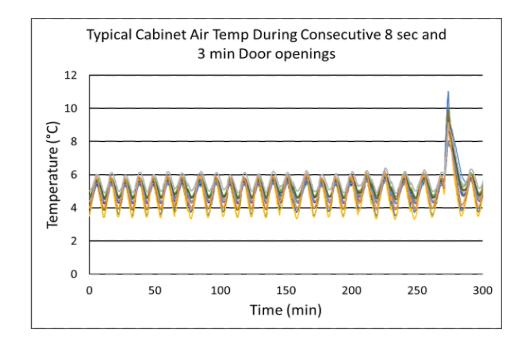

Controller, Configuration, Alarms and Monitoring				
Controller technology	Parametric, microprocessor, LED display with 0.1°C resolution			
Temperature setpoint range	1°C to 10°C (Setpoint must remain unaltered from the factory setting to remain compliant with NSF/ANSI 456 Standard for Vaccine Storage requirements)			
Display probe	Calibrated, stainless steel			
External alarm connection	State switching remote alarm contacts			
	Visual and audible indicators			
Alarms	High / Low temperature, compliant with alarm requirements defined in the NSF/ANSI 456 Standard for Vaccine Storage			
Simulator ballast	Glass bead thermal media			

Performance data acquired at 22°C ambient, using NSF/ANSI 456 compliant validation ballast probes, empty chamber, during stabilized steady state operation and a DAQ sampling rate of one measurement every 10 seconds

- 1 Uniformity is defined as the maximum variance in temperature across all probes at any point in time over the testing period
- 2 Stability is defined as the maximum variance in temperature experienced by any single probe over the testing period

3 - Temperature performance for all loaded and unloaded door opening protocols, all alarm, controller and probe requirements as defined in the NSF/ANSI 456 standard for vaccine storage

4 - Data per Energy Star test results or equivalent testing and calculation. Heat rejection based on daily averages, not continuous operation. Performance exceeds Energy Star requirements.




Images

Product Data Sheet

Undercounter 4.6 cu. ft. Built-in Glass Door Vaccine Refrigerator ADA - Certified to NSF/ANSI 456 Standard for Vaccine Storage



Dimensions					
	Width	Depth	Height	Door Swing	Total open Depth
Exterior	23 7/8"	26"	31 15/16"	23 1/2"	46"
Interior	19 1/4"	17 1/2"	22"		

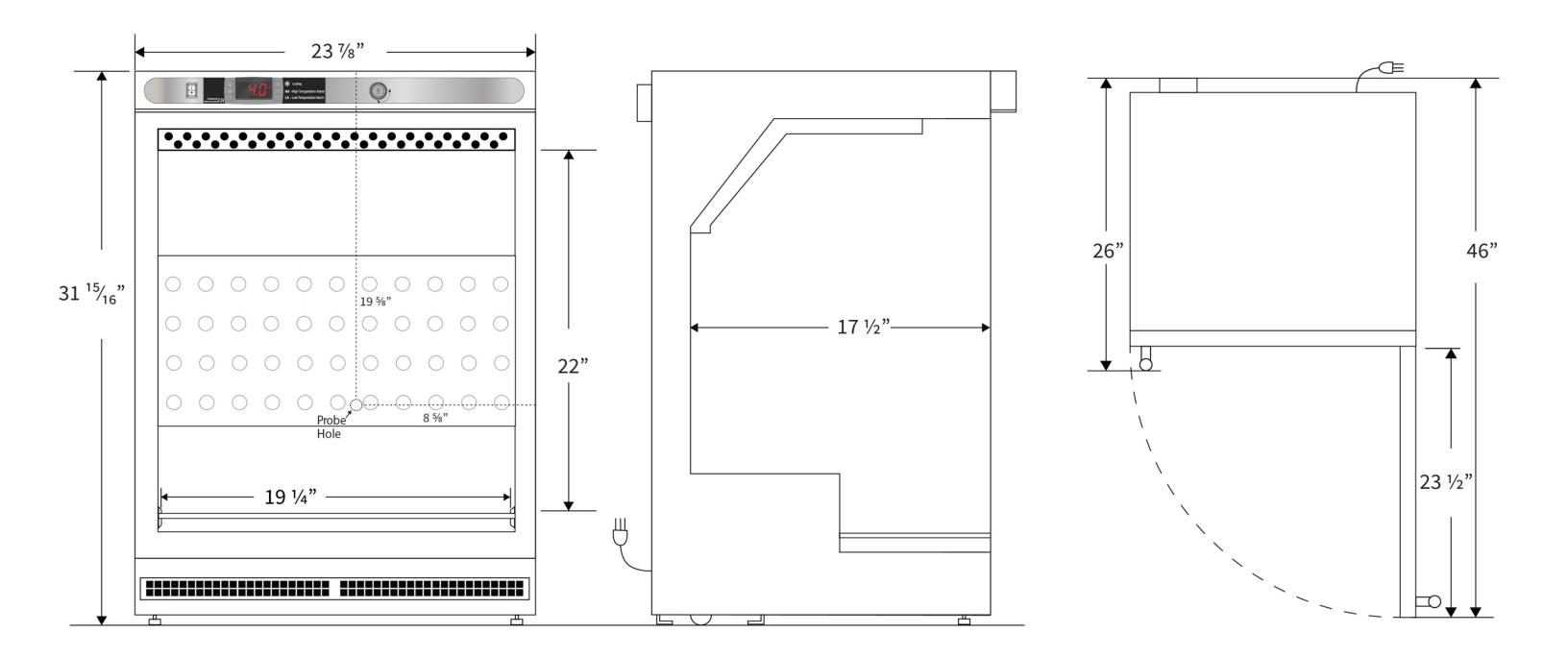

Contact		
Customer Service	800-648-4041 Option 3	customerservice@horizonscientific.com
Technical Service	800-648-4041 Option 5, Parts Option 4	technicalservice@horizonscientific.com
Rev_09302022		